

MEM,NEAA(With HEPES) Complete Medium

Product manual

Basic Information

Cat.NO	Size	Shelf	Form	Storage	Transportation
CMC0022	500mL	6 months	Liquid	Store at 2-8°C away from light	2-8°C

Product Introduction

Components	MEM,NEAA(With HEPES)(CMB0063)+10% Nutrient+1% Supplement
Concentration	Ready-to-use
Bacteria Detection	Negative
Fungi Detection	Negative
Mycoplasma Detection	Negative
Endotoxin Content (EU/mL)	< 3

Background

MEM medium (Minimum Essential Medium) is also called the minimum essential medium, minimum basic medium or low-limit Eagle medium. It was developed by Harry Eagle based on Eagle Basic Medium (BEM). It is the most basic and widely applicable medium and one of the most commonly used mediums in animal cell culture. MEM medium contains only 12 essential amino acids, glutamine and 8 vitamins. It has simple ingredients and is mainly used for the culture of adherent cells. It can also be used for other types of cell culture after the formula is modified. MEM medium containing NEAA (non-essential amino acids) is based on MEM medium and adds 7 NEAAs: L-alanine, L-glutamic acid, L-asparagine, L-aspartic acid, L-proline, L-serine and glycine. It can reduce the side effects of the production of non-essential amino acids by cells themselves during cell culture and effectively promote cell proliferation and metabolism.

HEPES is an excellent biological buffer with no toxic effect on cells. The culture medium with HEPES added can maintain a constant pH range for a long time, which can effectively prevent the adverse effects of large pH fluctuations in the culture medium on cell growth.

The culture medium with NEAA (non-essential amino acids) is based on the original culture medium and added with seven NEAAs: L-alanine, L-glutamic acid, L-asparagine, L-aspartic acid, L-proline, L-serine and glycine. It can reduce the side effects of the cells' own production of non-essential amino acids during cell culture and effectively promote cell proliferation and metabolism.

Precautions

1. During the entire process, be sure to pay attention to aseptic operation to avoid contamination;
2. To maintain the best use effect of this product, please do not perform freeze-thaw treatment;

-
3. This product is only used for research or further research, not for diagnosis and treatment.
 4. It is not suitable for long time storage at room temperature.
 5. This product is a ready-to-use medium. If there is no special need, don't add serum, penicillin and streptomycin. It can be used directly.